

Colifast Field Kit: Presence/Absence Test

Rough Guide


1. Extract water sample




2. Add to prefilled bottle




3. Incubate for 12 hours at 37 °C

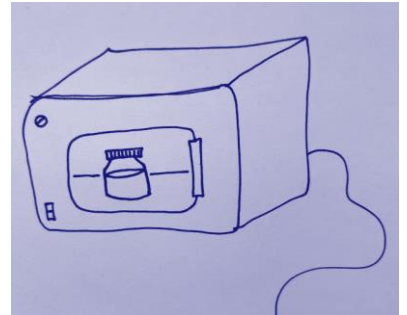


4. Subsample to cuvette

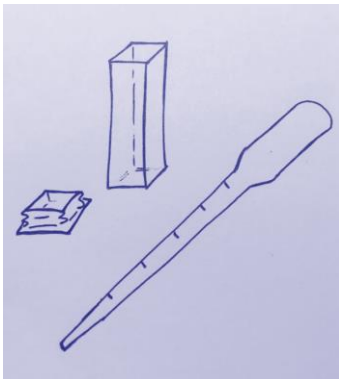

5. Add 3 drops of sample developer

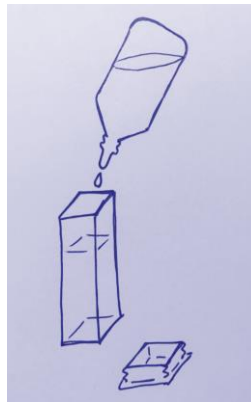

6. Analyse subsample with Colifast fluorometer

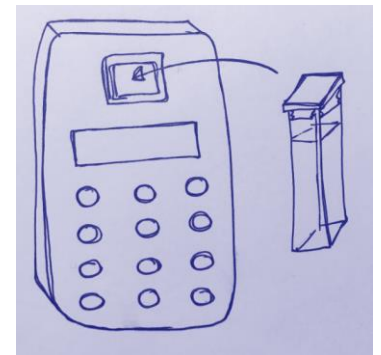

7. Value above 200 indicates presence of total coliforms (> 1 TC / 100 mL)

COLIFAST®

Technical support: Colifast AS, Strandveien 33, P.O. Box 31, 1324 Lysaker, Norway.

Tel: +47 67100510, Fax: +47 67100520, e-mail: post@colifast.no, web-site: www.colifast.com