


Colifast Field Kit: Rapid Test Rough Guide


1. Extract water sample






2. Add to prefilled vial

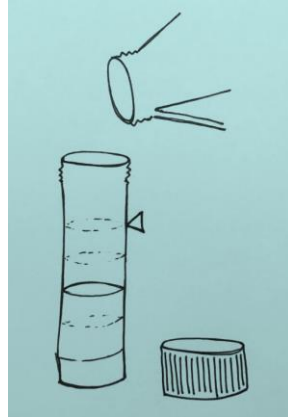

3. Incubate at 37 °C

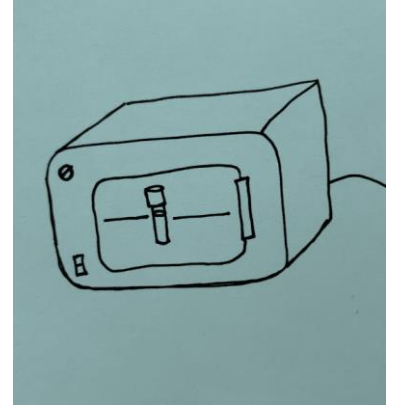

4. After 15, 45 and 75 min: Subsample and analyse with NaOH addition using Colifast Microdetector

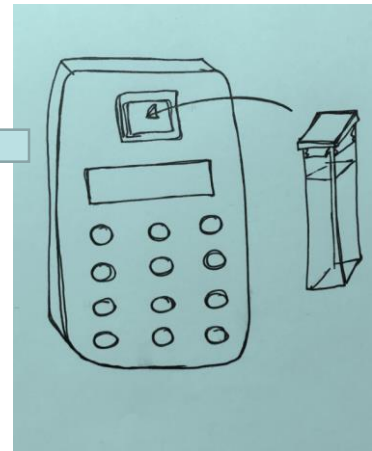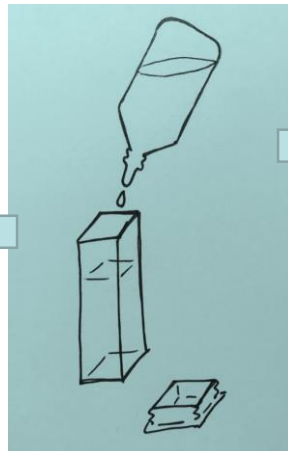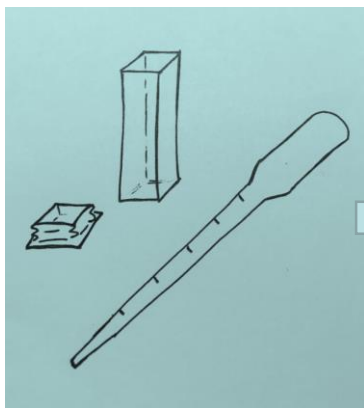

CMD slope (MU Production/hour)	cfu/100 mL
<2,5	~ 0
2,5	500
5	1 000
25	5 000
50	10 000
100	20 000
500	100 000

1 slope unit = 200 cfu / 100 mL

5. Calculate slope

COLIFAST[®]

Technical support: Colifast AS, Strandveien 33, 1366 Lysaker, Norway.

Tel: +47 67100510, Fax: +47 67100520, e-mail: post@colifast.no, web-site: www.colifast.com