


Colifast Field Kit: Quantitative Test (MPN) Rough Guide


1. Extract water sample




2. Add to pre-filled vial

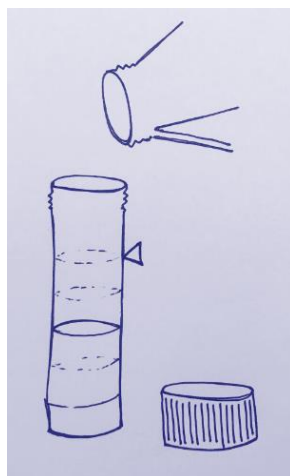


3. Subsample to cuvettes

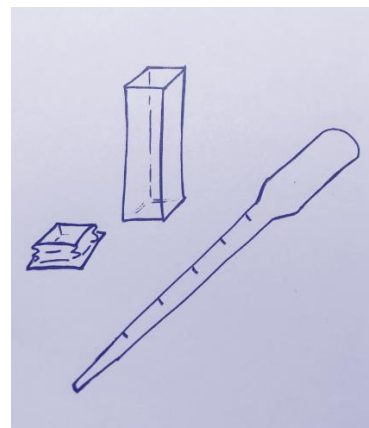


4. Incubate for 11 hours
at 37 °C

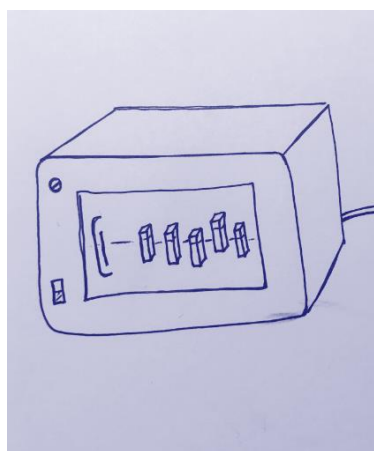

5. Add 3 drops of
sample developer

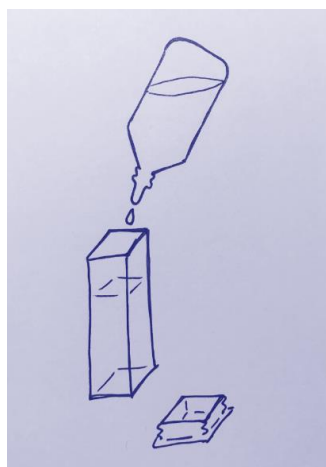

6. Analyse subsamples
with Colifast fluorometer

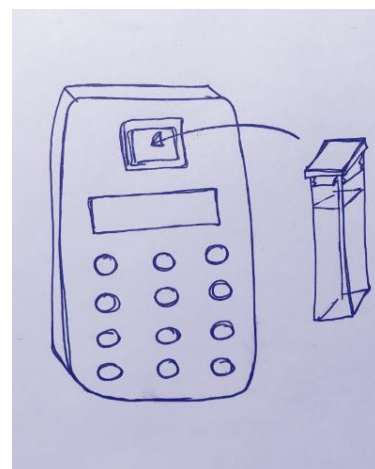

Number positive cuvettes (of 6)	# <i>E. coli</i> pr 100 ml
0	UNDER
1	12
2	27
3	47
4	77
5	136
6	OVER

7. Use table to determine
the Most Probable Number

COLIFAST®

Technical support: Colifast AS, Strandveien 33, P.O. Box 31, 1324 Lysaker, Norway.

Tel: +47 67100510, Fax: +47 67100520, e-mail: post@colifast.no, web-site: www.colifast.com